


# LIEBHERR

## Scientific and Healthcare Refrigerator and Freezer Quick Guide

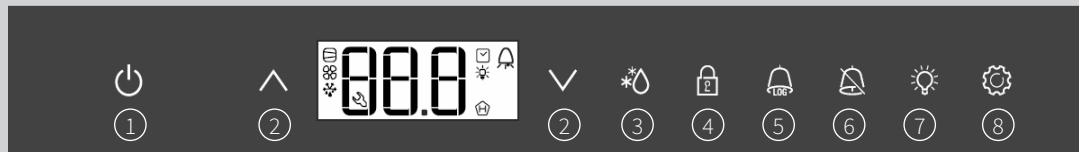

- |   |  |
|---|--|
| ① ON/OFF button                                   | ② Selection buttons                            |
| ③ Defrost button (for manual defrost)             | ④ Keypad lock                                  |
| ⑤ Button for calling up stored alarm events (log) | ⑥ Audible alarm OFF button & for exiting menus |
| ⑦ Interior light ON/OFF button                    | ⑧ Enter/Set button                             |

### Check temperature setpoint

5°C (for vaccine or medicine storage)

Press and hold . Temperature display flashes.

Adjust number to desired number by using or .

Press the to save the setting.

### Check temperature alarm

3 (for vaccine or medicine storage)

Press and hold down . Display = **r5**.

Press the or . Display = **AH** (high alarm limit) or **AL** (Low alarm limit).

Press . Number shows on display.

Adjust the number to desired temperature difference by using or .

AL or AH is designed to be set to temperature difference between temperature setpoint (5°C for vaccine or medicine storage) and high/low alarm temperature (8/2°C for vaccine or medicine storage). Hence, 3 for vaccine or medicine storage.

Press the to save the setting.

### Read high and low temperature log

Press and hold down . Display = **HAH**.

Press Display = **rL**.

Press . Read low temperature in °C since the last read and reset.

Press to return to **rL**.

Press . Display = **rH**.

Press . Read high temperature in °C since the last read and reset.

Press to return to **rH**.

### Read high and low temperature log

Press and hold down . Display = **HAH**.

Press until the display shows **rL**.

Press . Display = **0-999**.

Press and hold down . Display = **rE5**. Display = **0**.

### Temperature and power failure alarms

flashes & HF on display

- A power failure occurred.

flashes & HA & HI on display

- A high temperature alarm occurred.

flashes & HA & LO on display

- A low temperature alarm occurred.

### Clear temperature (power failure) alarms

Press and hold down . Display = **HAH**.

(Press until **HFH** appears on the display to clear power failure alarm).

Press and hold down and at the same time, Display = **rE5**.

### Read temperature alarm log

Press and hold down . Release the button until display shows **HAH**.

Press . Display = **1** (1 temperature alarm has occurred).

Press . Return to **HAH**.

Press . Display = **HA** Last temperature alarm

Press . Display = **900** (year)

Press . Display = **100** (month 1-12)

Press . Display = **100** (day 1-31)

Press . Display = **h00** (hour 0-23)

Press . Display = **m00** (minute 0-59)

Press . Display = **t00** (alarm duration in minutes)

**Note: To exit any setting and return to normal display, press and hold down for approximately 5 seconds.**

For further assistance please call Andi-Co Australia on 03 9569 1255 or email [service@andico.com.au](mailto:service@andico.com.au)

Andi-Co Australia

1 Stamford Road, Oakleigh, VIC 3166  
[andico.com.au](http://andico.com.au) | [home.liebherr.com.au](http://home.liebherr.com.au)